


脂肪酸甲酯的顺反异构

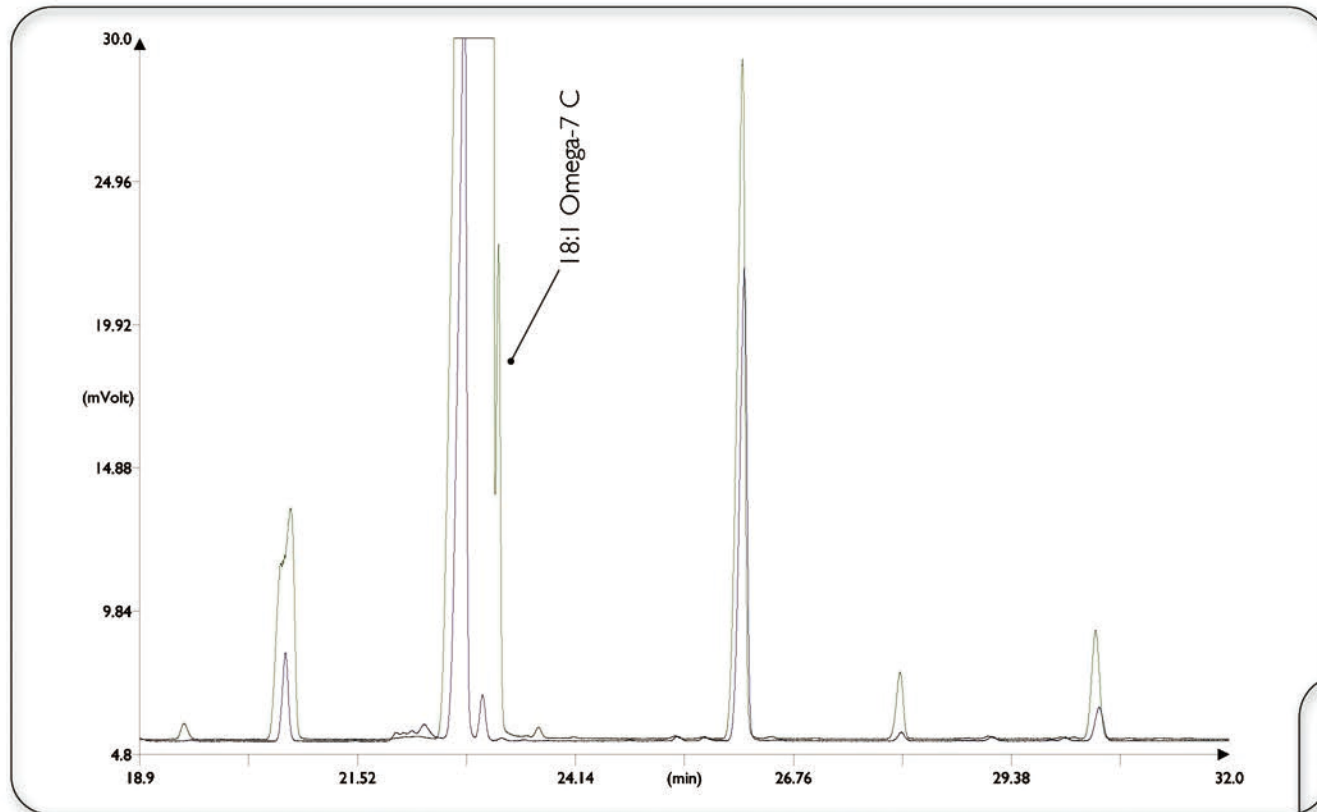

Column: **MEGA-10 - 0.25mm, 0.20 μ m, 60m**
Catalog Code: C-10-025-020-60

条件:

进样口: 分流 230 $^{\circ}$ C, 35mL/min分流比, 0.5 μ L进样量

检测器: FID 250 $^{\circ}$ C

柱温箱程序: 150 $^{\circ}$ C, 1 $^{\circ}$ C/min

载气: 氦气, 150kPa

Sample:

— MEGA-10 Test Mixture: calibrated mix of *cis/trans* FAMES.

— Real sample overlapping.

